

# GENERAL USE INSTRUCTIONS

## FS SERIES

### PRODUCT REFERENCE

FS series

### PRECAUTIONS

- Do not open the FS Series devices or FLOWBOARD
- Use only the cables cables provided by Fluigent
- Prevent foreign objects from entering the FS Series and FLOWBOARD
- Place the product in a stable location with a level surface and a stable support
- Respect the temperature compatibility (from 5°C to 50 °C)
- Filter your solutions. If possible, add a filter in the fluidic path and clean the flow sensor after each use, especially the XS models. The diameter of the XS capillary is small: 25 µm. Fluigent is not responsible in the event of clogging or surface modifications.
- Do not allow the flow sensors to dry with media in the flow path without flushing first.
- Fluigent advises to clean the sensors after use.
- The FS series yellow plugs must be connected for storage
- Verify the fluid compatibility with the FS Series wetted materials before using it or ask Fluigent customer support.

### SYMBOLS



**Disposal** Do not dispose of electric devices together with household waste material! In observance of European Directive 2012/19/EU on waste electrical and electronic equipment and its implementation in accordance with national law, electronic items that have reached the end of their life must be collected separately and returned to an environmentally compatible recycling facility.